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Originalversion

OPERATING MANUAL

MODE D'EMPLOI

MANUAL DE INSTRUCCIONES

D 82000 / 82100 / 82200

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WATER DISTILLING APPARATUS

APPAREIL A DISTILLER L'EAU

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Important notes:

The operating manual is part of the device. Before initial operation of the unit, please carefully read and observe the operating instructions and keep it. For safety reasons the unit may only be used for the purposes described in these present operating instructions. Please also observe the operating instructions for the units to be connected

All specifications in this operating manual are guidance values which are valid at the time of printing. However, for technical or commercial reasons or in the necessity to comply with the statutory stipulations of various countries, the manufacturer may perform additions to the unit without changing the described properties.

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Höhe 750 mm

Gewicht 11 kg

1	WARNING AND SAFETY INSTRUCTIONS	13
2	CONSTRUCTION OF THE DEVICE	14
3	PUTTING INTO OPERATION	15
4	OPERATION	16
5	CLEANING	16
6	PARTS.....	17
7	TECHNICAL DATA.....	18

Notes to the Manual

The provided manual will allow you the proper and safe handling of the product. For maximum security, observe the safety and warning instructions in the Instructions.

The pictogram  has the following meaning:

- Warning of a general danger to personnel and equipment.
- Non-compliance results (can result) in injury or material damage.

 Important information for device use.

 Refers to another part of the operating manual.

Status at time of printing

Advanced technology and the high quality of our products are guaranteed by a continuous development. This may result in differences between this operating manual and your product. We cannot exclude mistakes. We are sure you understand that no legal claims can be derived from the information, illustrations and descriptions.

A potentially more recent version of this manual is available on our internet website at www.si-analytics.com. The German version is the original version and binding in all specifications.

Guarantee

We provide guarantee for the device described for two years from the date of purchase. This guarantee covers manufacturing faults being discovered within the mentioned period of two years. Claim under guarantee covers only the restoration of functionality, not any further claim for damages or financial loss. Improper handling/use or illegitimate opening of the device results in loss of the guarantee rights. The guarantee does not cover wear parts, as lobes, cylinders, valves and pipes including the thread connections and the titration tips. The breach of glass parts is also excluded. To ascertain the guarantee liability, please return the instrument (see backside of this manual) and proof of purchase together with the date of purchase freight paid or prepaid.

Copyright

© 2016, Xylem Analytics Germany GmbH

Reprinting - even as excerpts - is only allowed with the explicit written authorization.

Printed in Germany.

1 Warning and safety instructions

This manual contains basic instructions which must be observed during installation, operation and maintenance of this device.

The Water Distilling Apparatus meets the requirements of Protective Class I.

It has been built and tested in accordance with DIN VDE 0700, Part 1/04.88 (HD 251 S3 and Changes 1-3) and DIN VDE 0700, Part 6/11.89 (HD 275 S1, Change 1) in connection with the draft of DIN VDE 0700, Part 6A8/04.89, Protective Measures for Electric Heating Appliances. It has left the factory in perfect technical and safety-related condition. In order to preserve this condition and ensure operation without danger, the user must observe the notices and warnings which are contained in these operating instructions.

 For safety and functional reasons, the Water Distilling Apparatus may be opened only by authorised persons; e. g. work at the electric appliance must only be carried out by trained specialists. With unauthorised operation of the device as well as careless or deliberate damaging, the warranty is no longer void.

 Before switching on, it is to be ensured that the operating voltage of the Water Distilling Apparatus and the network voltage are in agreement. The network plug connector may only be introduced into a socket with a protective contact. The protective effect may not be nullified through the use of an extension cable without a protective conductor. Any interruption of the protective conductor within or external to the Water Distilling Apparatus or the loosening of the protective conductor connection can lead to a situation in which the Water Distilling Apparatus represents a danger. Intentional interruption is not allowed.

 If it has to be assumed that safe operation is impossible, the Water Distilling Apparatus has to be put out of operation and secured against inadvertent putting to operation. (In this case please switch the unit off, pull plug of the mains cable out of the mains socket, and remove the unit from the place of work).

Examples for the assumption that a safe operation is no longer possible,

- the package is damaged,
- the unit shows visible damages,
- the unit does not function properly
- liquid has penetrated into the casing
- if the unit has been altered technologically or if unauthorized personnel tried or succeeded to open the instrument as attempt to repair it.

 Water Distilling Apparatus whose current carrying capacity is greater than 16 amps, are supplied without wall plugs. Connection to the mains supply is to be carried out by an authorised electrician, and the power consumption indicated on the identification plate should be observed.

 The Water Distilling Apparatus must not be stored or operated in humid rooms.

2 Construction of the device

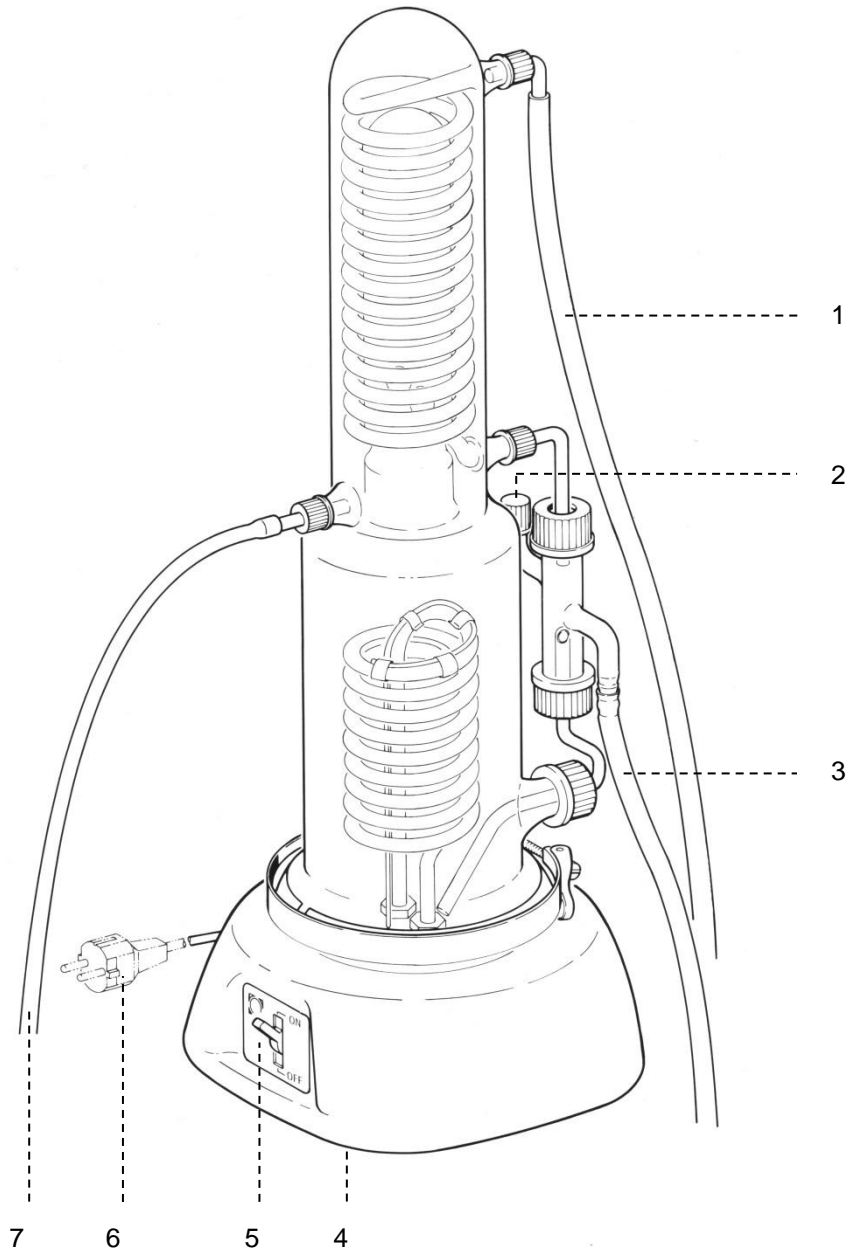
Fig. 1

- 1 Water inlet tube (to be connected to water main)
- 2 Venting cap (to be removed before putting into operation)
- 3 Water outlet
- 4 Reset button in base plate (overheating protection)
- 5 Master switch
- 6 Mains supply (observe connected load values)
- 7 Distillate

3 Putting into operation

1. Connect Water Distilling Apparatus as shown in Fig. 1
2. Adjust regulating part DZ 8103 according Fig. 2

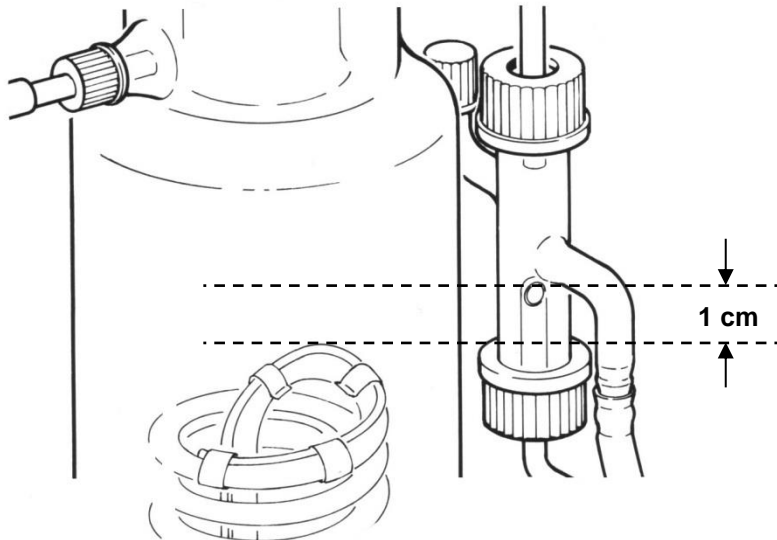
Fig. 2

i The marking lines drawn show the correct adjustment of the DZ 8103 control component, required for water level control in the water distillation apparatus:

The water outlet lower edge is required to be located at a level **1 cm** higher than the heater coil upper edge! (This ensures the heater coil is completely immersed in water)

i In the event of the water level falling to below the heater coil upper edge, the apparatus is disconnected!

3. Remove venting cap (2). Check for tight fit all other plastic screw caps and, if necessary, re-tighten.
4. Open tap water valve (water faucet):

! Before switching on Water Distilling Apparatus, fill boiling vessel with water

5. As soon as water flows from water outlet (3) (after approx. 2 minutes) turn on heating with the master switch (5).
6. Distillation will start after approx. 5 minutes.

4 Operation

The Water Distilling Apparatus will have attained optimum working conditions when temperature of out flowing water (3) is of approx. 60 °C. Distillate (7) running off is then free from CO₂ and loss of energy is small.

The Water Distilling Apparatus is provided with overheating protection:

 In case of water shortage in boiling vessel, heating turns off automatically. After restoring normal working conditions, distiller will again be ready for use after **pressing hard** on reset button (4).

 Press reset button of overheating protection **only after** Water Distilling Apparatus has **cooled down!**

To obtain uniformly good water quality, it is necessary to allow the salt-enriched water in the boiling vessel to run off daily. To this effect, secure again venting cap (2), after having turned off the heating, open water faucet a little more for a few seconds and then shut off completely. Boiling vessel content will siphon off.


5 Cleaning

 Cleaning with a detergent solution is necessary after 1 to 4 days of operation according to the salt content of the tap water:

1. Drain Water Distilling Apparatus.
2. Remove venting cap (2) and insert funnel
3. Slowly fill in 0.5 l detergent solution:
avoid heavy foaming, if necessary, interrupt filling operation
(1 l detergent solution Type No. DZ 8050, 5 l detergent solution Type No. DZ 8053)
4. Re-place in position venting cap (2) and tighten again.
5. Close off drain hose (3) (with pinch cock, for instance)
6. Refill boiling vessel with tap water up to approx. 1 cm above heating coil.
7. Heat up to boiling point temporarily, avoid heavy foaming.
8. Wait until water incrustation (boiler scale) has dissolved up to filling level.
9. Fill boiling vessel completely with tap water, so that the last boiler scale residue may also dissolve.
10. Open drain hose (3). Boiling vessel content will siphon off.
11. Close off drain hose (3). Fill again boiling vessel completely with tap water and allow to siphon off.
 We recommend to **repeat** the washing process **five times**.
12. Do not use distillate after putting the Water Distilling Apparatus into operation, until the conductivity or the pH-value of the distillate have attained the value required. Without measuring, we recommend not to use the distillate of the first hour of operation.

 The Water Distilling Apparatus may never be submerged into water!

6 Parts

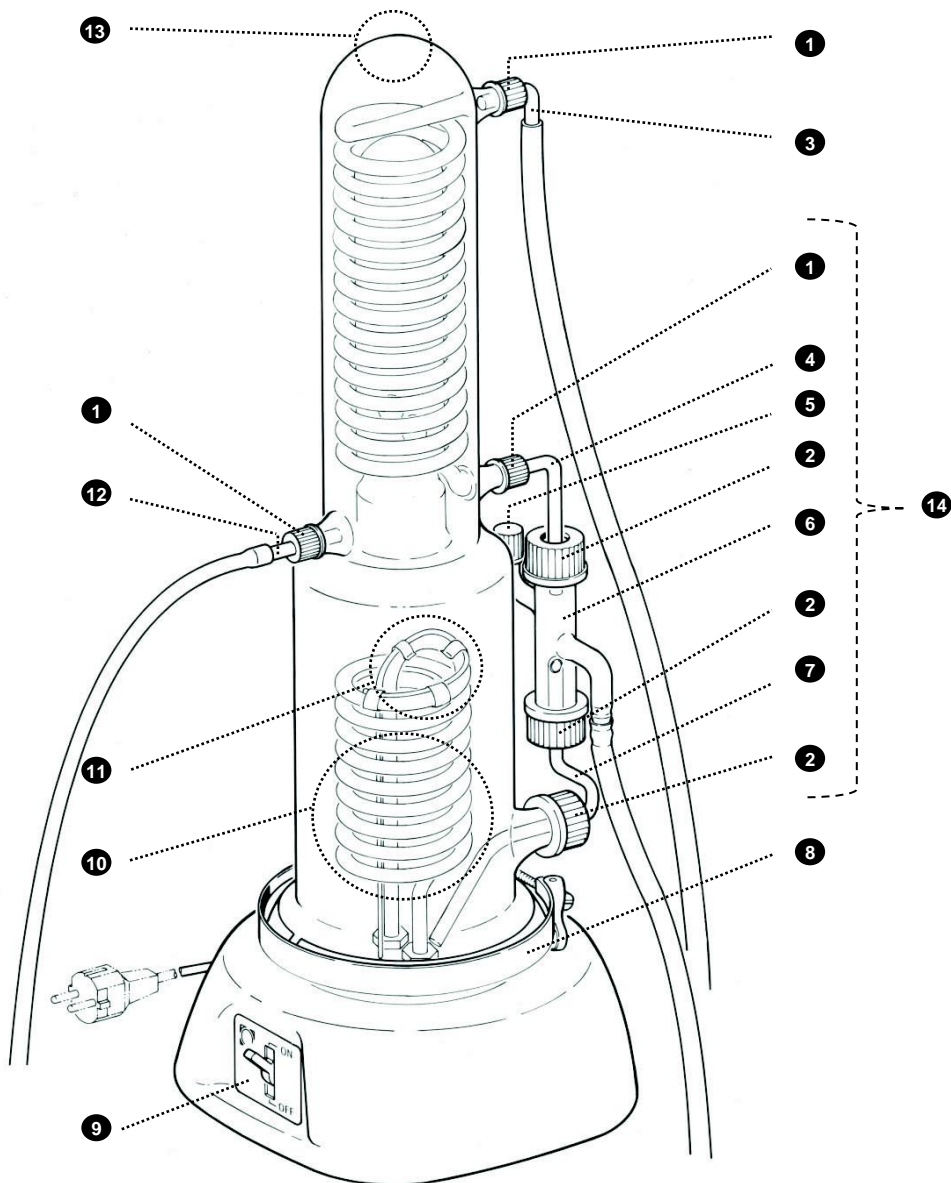1 DZ 8215 (cap) +
DZ 8206 (seal)

2 DZ 8208 (cap) +
DZ 8218 (seal)

3 DZ 8104

4 DZ 8102

5 DZ 8411

6 DZ 8103

7 DZ 8106

8 DZ 8757 + DZ 8257

9 DZ 8524 (for D 82000)
DZ 8505 (for D 82100/82200)

10 DZ 8500 (4I/230V for D 82000)
DZ 8600 (7I/230V for D 82100)
DZ 8700 (7I/380V for D 82200)

11 Overheat protection
DZ 8701
DZ 8702

12 DZ 8105

13 DZ 8101 (for D 81100/81200/
82000/82100/82000)

DZ 8107 (for D 82100/82200)

14 Level controller (complete)
composed of:

1x DZ 8102

1x DZ 8103

1x DZ 8106

1x DZ 8206

3x DZ 8208

1x DZ 8215

3x DZ 8218

7 Technical data

Type-No.	Electrical input power V/A	Output H₂O dist. l/h
D 82000	230/14	4
D 82100	230/23	7
D 82200	400/3x8	7

Type of heating	resistance heating
Dimension of base	300 x 300 mm
Height	750 mm
Weight	11 kg

**EU – KONFORMITÄTSERKLÄRUNG / EU – DECLARATION OF CONFORMITY
UE – DECLARATION DE CONFORMITE / UE – DECLARATION DE CONFIRMIDAD**

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Xylem Analytics Germany Sales GmbH & Co.KG